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YC-015 LAB SPRAY DRYER

YC-015实验型喷雾干燥机是上海雅程仪器设备有限公司与美国PILOTECH公司在实验喷雾干燥领域多年合作开发的结晶。整机设计紧凑，可安放在专门设计的不锈钢机架或实验台上，自成一體，无需其他设施即可运行。一键式开机，彩色液晶触摸屏操作，可采用完全自动或人工监控两种运行模式，方便操作和实验过程的监控。YC-015, made by Shanghai Yacheng Inc. cooperating with PILOTECH Inc on Spray drying research for many years, is self contained and supplied complete and ready for immediate operation. All major components are housed within a stainless steel cabinet and the unit can be used on a bench top or with an optional stainless steel stand. The system was designed to be stand by a keyboard, conducted by a colorful crystal screen of touch guidance mode, and allowed two modes of run: Automatic-mode, and Eye-monitored mode for the purpose of easily controlling experimental process.

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手动控制：可方便从自动状态切换至手动状态来修改参数
Two modes of run: Automatic-mode, and Eye-monitored mode for the purpose of easily controlling experimental process. PLC controller, LCD touch panel control fascia with flow path of system.

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With cooling device for temperature-sensitive material (optional)

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Shutdown protection: press the stop button and the machine (except the fan) stops working immediately so that heating parts will not be burnt due to misuse (forced shutdown of the fan).

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Built-in air oilless compressor, the noise is low than 60db

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Evaporation rate of water at inlet temperature of 250°C using Standard: Approximately 1500 ml/hour~ 2000ml/hour

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Spray chamber, cyclone separator, receiving tank are all made of Borosilicate glass which can resist strong acid/alkali, high temperature. It can work in a no-pollution and stable environment, and the whole process can be inspected. All the spare parts are easy to install and wash.

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As standard a touch screen to display temperature, airflow volume, air pressure, pump speed and de-blocker frequency.

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To assure accurate temperature control and easy parameter adjusting, YC-015 applies PID. Temp. Precision: ±1°C, dry cycle time: 1.0-1.5S

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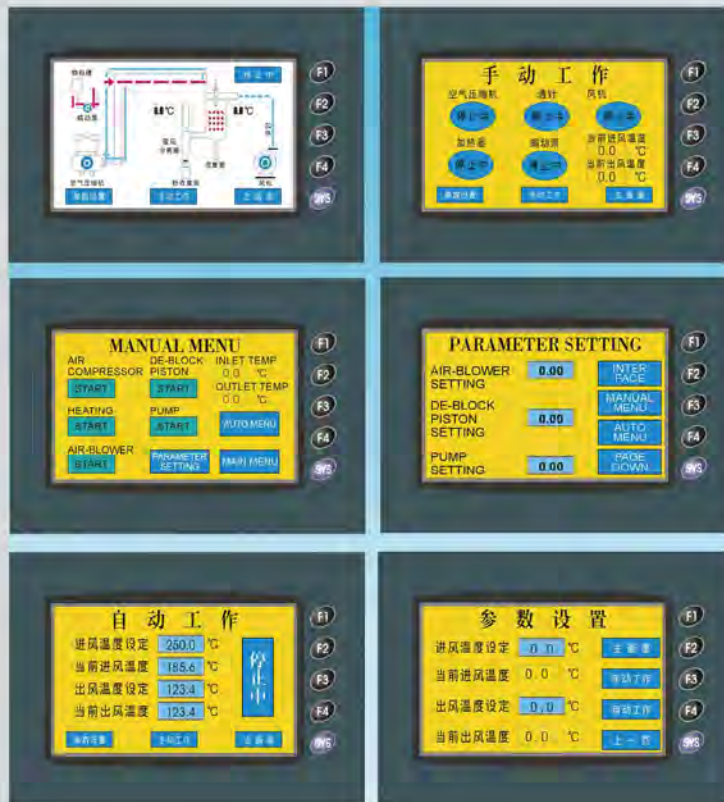
Two modes of run: Automatic-mode, and Eye-monitored mode for the purpose of easily controlling experimental process . PLC controller, LCD touch panel control fascia with flow path of system.

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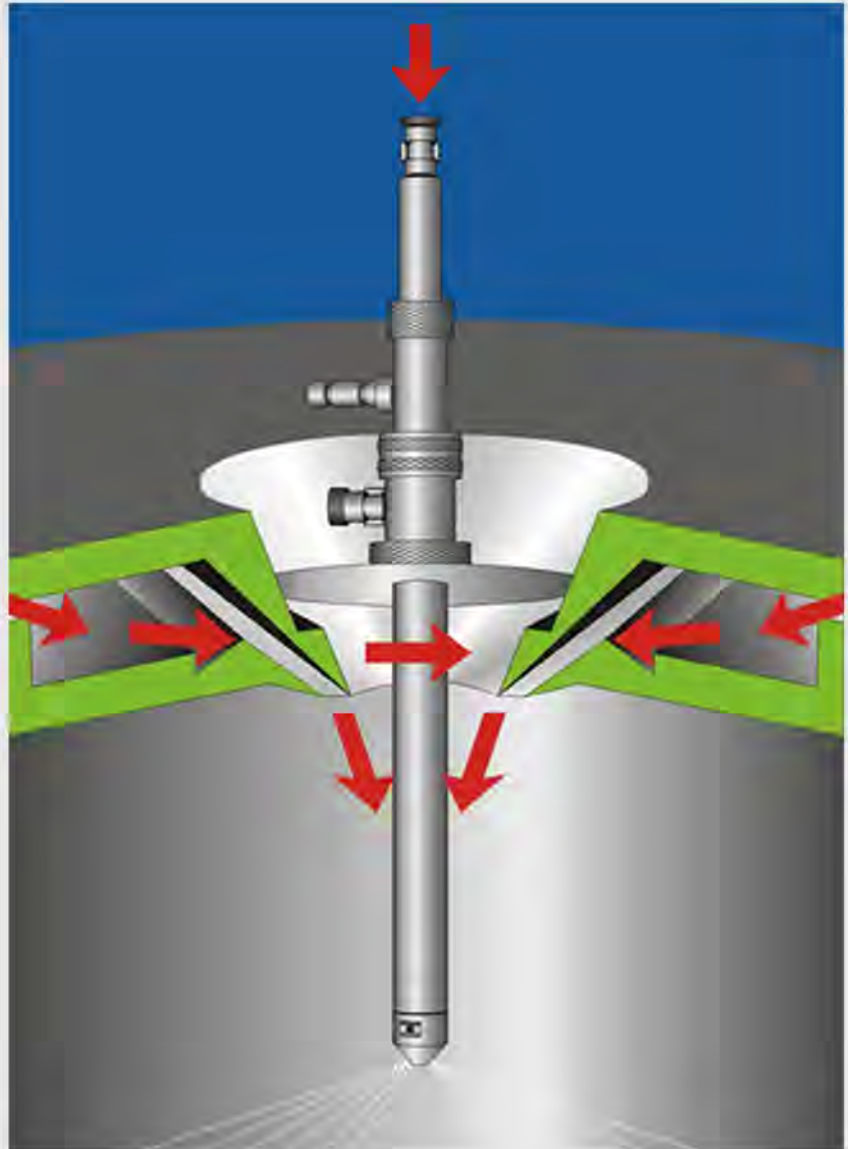
Spray chamber, cyclone separator, receiving tank are all made of Borosilicate glass which can resist strong acid/alk li, high temperature. It can work in a no-pollution and stable environment, and the whole process can be inspected. All the spare parts are easy to install and wash.

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A filter has been equipped to the inlet unit to prevent the sample from contamination.

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The self-priming peristaltic pump delivers the sample liquid from a container through a small diameter jet into the main chamber to avoid secondary pollution.. To make sure the process conducted steady and smoothly, magnetic agitator is optional to make the solution well mixed.

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Narrow size distribution of particles.

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Automatic de-blocking device prevents the nozzle from becoming blocked and variably controlled